

Disposable Sack Sampler

Take a sample through the side of a Sack

The disposable Sack Sampler is a hygienic, single use device that has been specially designed to take samples through the side of a sack.

- Individually Bagged
- Made from FDA, EC 1935/2004 & EU 10/2011 Conforming Materials
- BSE/TSE Free
- Full Batch Traceability
- Suitable for use in ATEX areas
- Available sterile (SAL 10⁻⁶) and non-sterile

Operation:

1. Insert the sampler through a sampling hole in the side of the sack.
2. Angle sampler downwards to allow the product to flow into the sample collecting bag
3. Seal hole with an Ultra Adhesive label



Part No.	Length	Diameter	Material	Sterile	Quantity per Box
8085A-100	400mm	25mm	HDPE	No	100
8085A-100S	400mm	25mm	HDPE	Yes	100
8085N-100	400mm	14mm	HDPE	No	100
8085N-100S	400mm	14mm	HDPE	Yes	100
8080A-450	Sample Collecting Bag 450ml - for hygienic use - sterile			Quantity: 1000 bags	
8080A-900	Sample Collecting Bag 900ml - for hygienic use - sterile			Quantity: 1000 bags	

SteriWare® Sack and Bag Sampler

Perfect for taking samples from sacks and bags

New

The SteriWare Sack and Bag Sampler is perfect for taking samples from sacks and bags. This sampler is ideal for fast quality checks through the sides of the sacks. The SteriWare Sack Sampler is available in two diameters. The narrow sack sampler (25mm) is ideal for free flowing powders and for small products like pepper, corns and small grains whereas the large diameter sampler (45mm) is more suited to sampling larger products like nuts and beans.

The SteriWare Sack Sampler has an ergonomic handle and specially designed blade.

During sampling the operator needs to hold a bag or bottle over the rear end of the sampler to catch the powder or granules as they slide through the sampler.



Operation:

1. Insert the sampler through a sampling hole in the side of the sack.
2. Angle sampler downwards to allow the product to flow into the sample collecting bag
3. Seal hole with an Ultra Adhesive label

Part No.	Sterile	Length mm	Diameter mm	Material	Box Quantity
8688H-330	No	325	33	HDPE	100
8688H-330S	Yes	325	33	HDPE	100
8688J-330	No	325	45	HDPE	100
8688J-330S	Yes	325	45	HDPE	100

Use Ultra Adhesive Labels to seal the sacks after sampling.



See the full range of labels at: www.sampling.com